

INSTRUCTION MANUAL

CAT. 60610 Zerostat Anti-Static Gun



The Zerostat Anti-Static Gun produces negative and positive ions that couple with the positive and negative static charges on an object's surface, effectively neutralizing them.

Instructions:

1. Hold the Zerostat Anti-Static Gun about 300mm (12 inches) away from the object.
2. Squeeze the trigger slowly; a powerful stream of positive ions projects over a spread of about 400mm (16 inches).
3. Release it slowly, and negative ions are produced.

The Zerostat Anti-Static Gun lasts for at least 50,000 trigger operations, and requires no power supply, batteries, or refills.

Electron Microscopy Sciences

P.O. Box 550

1560 Industry Road Hatfield, PA 19440

TOLL FREE: 1-800-523-5874
TEL: 215-412-8400 WEB: www.emsdiasum.com
FAX: 215-412-8450 EMAIL: sgkcck@aol.com